



How long will silanized vials last on the shelf prior to use – FAQ

Sealed packs of silanized vials can be stored indefinitely without any practical limitations. However, once the packaging is opened, the vials' surfaces become susceptible to particle adsorption.

This is particularly true for particles from cardboard packaging or the surrounding air. The silanized surface of the vials tends to attract and hold these particles, which can render the vials unusable.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 09-27-2024