

Can I backflush the Cogent column filter for longer in-service life - FAQ

The Cogent column filter can be backflushed but...

Backflushing might push the internal filter element out of place. And not having a clog in the frit, but with slightly dislodged filter element suggests that filter is not working. Therefore, backflushing is not supported.

Click [HERE](#) for ordering information and a picture of the Cogent column filters.



Printed from the Chrom Resource Center

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 07-22-2024