

Rescaling graphs in the MicroSolv HPLC PQ Kit Excel template – Tips & Suggestions

Steps to manually re-scale the Y-axis in the PQ Kit template

Note, Residual graphs are not automatically set to scale of the data entered. As such, the Y-axis default settings may be out of ratio for the data values entered. By manually re-scaling the Y-axis, it will reflect your values for a more effective visual ratio.

Please use the following instructions to format and scale the Y-axis of graphs in the MicroSolv HPLC PQ Kit Excel template :


1. In the detector linearity tab of the PQ RID spreadsheet, right click on the detector linearity residual graph.




2. This will open a tab of options. Click **"Format Plot Area."**



3. This opens a new grey tab to the right of the screen, labeled, **"Format Plot Area."** Click the blue arrow next to the **"Plot Area Options."**

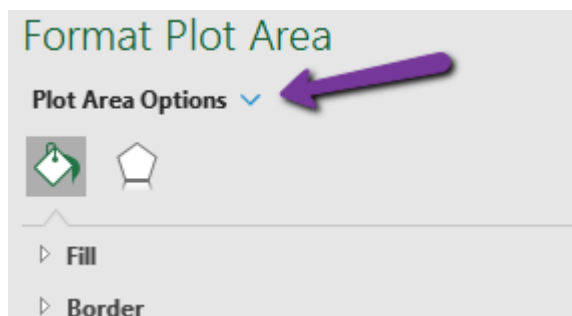

4. Next, click the 'Graph' icon.

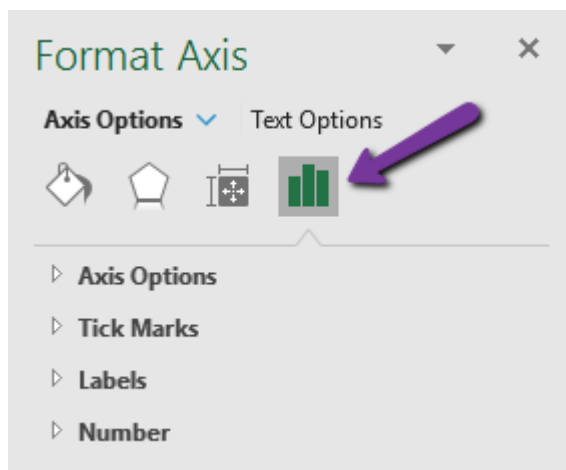

5. This will open a list under **"Plot Area Options."** click the **"Vertical (Value) Axis"** option.



6. This will open a tab to “**Axis Options.**” Then Check “**Reset**” on both the *Minimum* and *Maximum* **Bounds**.



7. Then check “**Reset**” on both the *Major* and *Minor* **Units**.



8. Your new graph should now be to scale and display all data points.



MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com