



## Standard silicone pre slit sealing mat compared to solvent resistant mat - FAQ

### **Comparison of 96 well sealing mats, silicone rubber v. chemical resistant U-2D mats.**

Silicone rubber mats should only be used with pure aqueous and buffer solutions.

The solvent resistant mat can be used with methanol, acetonitrile and water / buffers. These should not be used with non-polar solvents or aggressive acid or base such as ethyl acetate.

[U-2D Silicone Mats Product Page](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)