



## What are the temperature and pressure limits for amber glass autosampler vials - FAQ

**Amber glass vials are manufactured with 51-expansion, borosilicate glass, and according to the recommended exposure this glass is rated to 230°C for normal service or 490°C for extreme service.**

Please always verify the temperature limits of caps and septa, as each material will have different specifications.

We do not recommend pressurizing autosampler vials above 1 psi, as they are designed to be used as sample containment for HPLC or GC applications and we do not recommend using them outside the range of normal suggested use.

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)