





Remdesivir and GS-441524

A mixture of polar and non-polar drug compounds in ANP

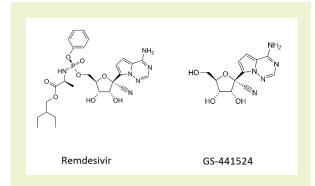

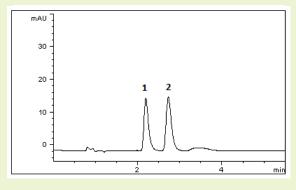

Notes: Remdesivir is a broad-spectrum antiviral drug that is able to diffuse into cells where it is converted to GS-441524 mono-phosphate. This achieved via the actions of esterases and a phosphoamidase; which in turn is further phosphorylated to its active metabolite triphosphate by nucleoside-phosphate kinases GS-441524. GS-441524 has previously been researched as a treatment for feline infectious peritonitis (FIP), a lethal coronavirus disease which affects domestic cats.



Column: Cogent Diamond Hydride™, 4µm, 100Å

Catalog No.: 70000-75P

Dimensions: 4.6 x 75 mm

Mobile Phase: A: DI H_2O + 0.1% formic acid (v/v) B: Acetonitrile + 0.1% formic acid (v/v)

Isocratic: 92% B

Injection vol.: 2µL

Flow rate: 1.0 mL/min

Detection: 254 nm

Sample: Stock SoluRemdesivir and GS-441524 in 0.1 mg/mL concentration

Peaks: 1. Remdesivir 2. GS-441524

Discussion

Remdesivir, an antiviral drug, and GS-441524, a pro-drug converted in the body as a ribonucleotide analog, is as a mixture that may be difficult to analyze in routine chromatographic analysis. As GS-441524 is polar, is makes it more difficult to retain in typical reversed phase conditions. By use of the Cogent Diamond Hydride[™] column in ANP, we have illustrated how this contrasting mixture of compounds can be retained and still possess excellent peak shape. This method demonstrates an easy HPLC method that can be easily transferred to LCMS.



9158 Industrial Blvd NE Leland, NC 28451 p: 1.732.380.8900 f: 1.910.769.9435 customers@mtc-usa.com www.Cogent-HPLC.com

APP-A-386