





Normal Phase Made Easier

Drying Solvents in the Mobile Phase not Necessary

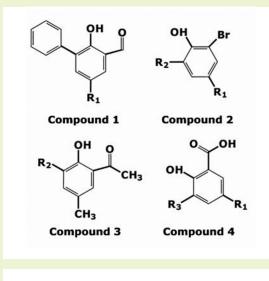

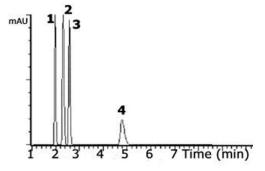

Notes:

1. This method can also be used with Atmospheric Pressure Chemical Ionization in the positive or negative mode: APCI+, APCI-

2. The column used for this application was previously used under Reversed Phase condition. Switching the column to Normal Phase mode required flushing 100% methanol for 10 minutes followed by 100% Methylene Chloride for 10 minutes. This process can be repeated over and over without damage to the column and can go back and forth between modes.



Column: Cogent Bidentate C18™, 4µm, 100Å

Catalog No.: 40018-75P

Dimensions: 4.6 x 75 mm

Mobile Phase: 95% Hexane/ 5% ethyl acetate

Injection vol.: 1µL

Flow rate: 1 mL/min

Detection: Photo Diode Array (UV)

Sample: 1 mg/mL of proprietary compounds were dissolved in the mobile phase

Peaks: 1. Phenolic compound with an aldehyde

- 2. Parent phenolic compound
- 3. Phenolic compound with a ketone
- 4. Phenolic compound with an acid group Photo Diode Array (UV)

Discussion

This Cogent Bidentate C18 column was used to quickly separate four proprietary phenolic compounds (precursors for a catalyst) under Normal Phase conditions. The separation shown is extremely reproducible (%RSD 0.2) and the solvents do not need to be dried before use as with other normal phase columns, saving time and lab resources.



9158 Industrial Blvd NE Leland, NC 28451 p: 1.732.380.8900 f: 1.910.769.9435 APP-A-06