



Plasma Extract

LC-MS method for polar compounds using UDC-Cholesterol



Method Conditions

Column: Cogent UDC-Cholesterol™, 4µm, 100Å

Catalog No.: 69069-05P-2

Dimensions: 2.1 x 50 mm

Solvents: A: DI H₂O / 0.1% formic acid (v/v) B: Acetonitrile / 0.1% formic acid (v/v)

adient:	time (min.)	%B
	0	90
	1	90
	7	20
	11	20
	12	90

Post time: 3 min

Gr

Injection vol.: 1 µL

Flow rate: 0.4 mL/min

Detection: ESI - POS - Agilent 6210 MSD TOF mass spectrometer

Sample: Proprietary lyophilized plasma sample was reconstituted in 60 microL 80% acetonitrile / 20% DI H₂O mixture

Peaks: 1. Urea 61.0396 m/z [M + H]+

- 2. Lactodifucotetraose 635.2393 m/z [M + H]+
- 3. Trimethylamine N-oxide 76.0757 m/z [M + H]+
- 4. L-carnitine 162.1125 m/z [M + H]+
- 5. Choline 105.1148 m/z [M + H]+

Discussion

When using the Cogent UDC-Cholesterol column and MS detection, it was possible to retain compounds present in a human plasma sample. The mechanism of retention of these polar compounds on Cogent UDC-Cholesterol column is unknown at this time. This Cogent column is known to retain and separate steroids and other compounds based on the shape of the molecules.

MICROS IV TECHNOLOGY

9158 Industrial Blvd NE Leland, NC 28451 APP-A-302