


## NMR structure confirmation data for Cogent UDC-cholesterol phases - Tech Information

The following solid state CP-MAS  $^{13}\text{C}$  NMR spectrum was recorded for a Cogent UDC-Cholesterol™ stationary phase in order to elucidate the structure. The assignments for the different carbon groups are shown below. The spectrum highlights the unique bonding of these phases, which is not possible with conventional silica based stationary phases.

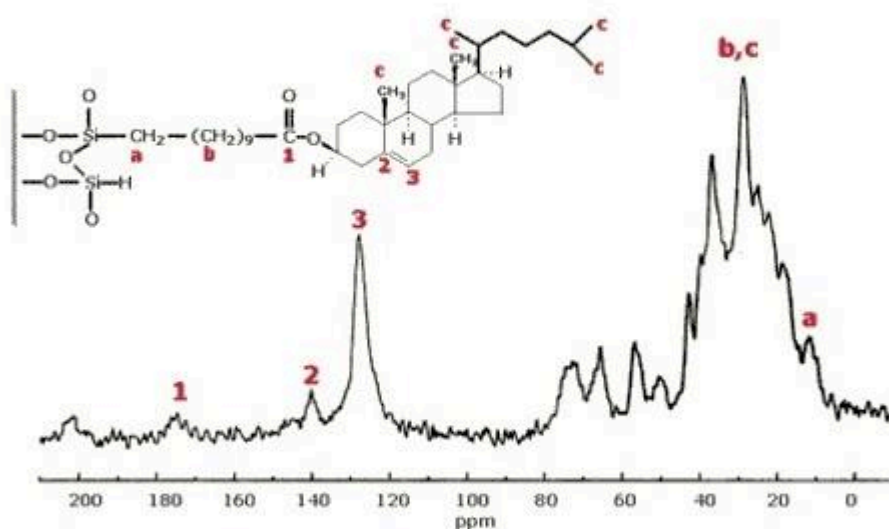

## UDC-Cholesterol

Click [HERE](#) for Cogent UDC-cholesterol HPLC column ordering information

Printed from the Chrom Resource Center  
Copyright 2025, All Rights Apply  
**MicroSolv Technology Corporation**  
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900  
Fax: (910) 769-9435  
Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)  
Website: [www.mtc-usa.com](http://www.mtc-usa.com)