



What is the purpose of using isopropanol IPA as part of the mobile phase solvent in ANP methods - FAQ

Some of our Cogent TYPE-C HPLC column application notes that feature ~~Aqueous Normal Phase~~ ANP use 50% isopropanol (IPA) as part of the aqueous component of the mobile phase. The main reason for this use is to help keep the column free of undesired adsorbed contaminants from a sample or sample matrix.

Contaminants can change the retention behavior of your ~~analyte~~ in any HPLC method and can also slowly elute off the column as carryover peaks in subsequent runs or as baseline noise and drift. We have found IPA (*or methanol can be used*) to be good at washing these contaminants off the column during the final part of an ANP gradient run. When performing HPLC or LCMS, adding IPA to the A (*aqueous*) solvent component will have little effect on the chromatography compared to using only aqueous-containing solvents, so it is something you should consider adding in the method before it has been fully developed.



Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900
Fax: (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com