

How can I get separation of tartaric and malic acid in a complex mixture? - FAQ.

Suppose you have a sample containing lactic acid, glycolic acid, formic acid, L-tartaric acid, DLmalic acid, and succinic acid. The Cogent Diamond Hydride™ HPLC column will be an excellent column choice for this separation. In the mobile phase that you will use, you will want to keep the acids ionized so that they will be more strongly retained on the column utilizing Aqueous Normal Phase (ANP) HPLC.

To do this, use a 10mM ammonium acetate <u>additive</u> in solvent A and 90/10 acetonitrile/10mM ammonium acetate in solvent B. For detection, use negative ion mode in LC-MS and look for the [M-H]- peaks in the EIC. For the gradient, start with the following:

Time (min) %B

0.0906.0309.03010.090

You may find that tartaric and malic acid co-elute under these conditions (peaks 4 and 5):



In this case, you can try a different gradient to get separation of peaks 4 and 5:

Time (min) %B

0.0	85
8.0	20
9.0	20
10.0	85



Click HERE for Cogent Diamond Hydride HPLC Column Ordering Information.



Printed from the Chrom Resource Center Copyright 2025, All Rights Apply **MicroSolv Technology Corporation** 9158 Industrial Blvd. NE, Leland, NC 28451 Tel: (732) 380-8900 Fax: (910) 769-9435 Email: customers@mtc-usa.com Website: www.mtc-usa.com