



## What is the best way to regain flow in the capillary if I allowed it to dry out overnight - FAQ

The best way to regain flow is to remove 1 or 2mm of each end of the capillary. Then gently force methanol through the capillary using a syringe pump.

Click [here](#) to view a convenient hand pump for this purpose.

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)