



What is the difference between RSA-Pro and the new RSA-Pro X surface treatments - FAQ

The biggest difference is that RSA-Pro™ surface treatment incorporated a vapor deposition type process that used ionic bonds with silanes to achieve hydrophobicity of the RSA glass and the RSA-Pro X™ surface treatment uses a very different, proprietary process of covalently bonding hydrophobic materials to the walls of RSA™ vials.

The RSA-Pro™ treatment is not very hydrolytically stable and could not be autoclaved or frozen and available only as a custom order item. RSA-Pro X™ will not lose its hydrophobic nature for years, even in the presence of water and they can be autoclaved and frozen and thawed effectively.



Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com