




## Can Snap Caps be used on Crimp Top Vials - FAQ

No, Snap Caps are specifically designed to be used only with vials that have a snap-ring finish—commonly referred to as [Snap Ring Vials](#). These vials feature a specially designed top that allows the "snap cap" to securely attach.

While crimp caps can technically be used on snap-ring vials, they are engineered to perform best with *serum finish vials*, also known as *crimp-top vials*. These vials provide the optimal sealing surface and neck structure for crimp caps, ensuring the best performance and seal integrity.

### Summary:

-  **Snap Caps** → Only for **Snap Top Vials** (with snap-ring finish)
-  **Crimp Caps** → Best used with **Crimp Top Vials** (serum finish)
-  **Do not use Snap Caps on Crimp Top Vials**

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)