



How do I store the Simplus brand bare silica CE capillaries - FAQ

For short term reuse, store Simplus™ capillaries at room temperature. Many users simply tape them to the wall near the instrument.

For long term storage and reuse, fill with water and store at room temperature. Before use, store at room temperature. Never refrigerate or heat these before use or during storage.

Click [HERE](#) for capillaries ordering information

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com