



## Can I use the MicroSolv Window Maker on all CE capillaries - FAQ

The MicroSolv Window Maker™ can be used on all polyimide coated, bare fused silica capillaries such as the Simplus™ or other brands.

It can be used on many internally coated capillaries but you should test the capillary to make sure that the coating is not removed by the high heat produced to remove the polyimide. If the coating is removed on the inside of the capillary, it could disrupt the flow and give erroneous results and therefore is not recommended. Contact the manufacturer of the CE capillary for more information.

Click [HERE](#) for MicroSolv Window Maker ordering information and pictures.

Printed from the Chrom Resource Center  
Copyright 2025, All Rights Apply  
**MicroSolv Technology Corporation**  
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900  
Fax: (910) 769-9435  
Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)  
Website: [www.mtc-usa.com](http://www.mtc-usa.com)