



## How do you replace the blade on the CleanCut90 PEEK tubing cutter tool - How To

There is a stainless steel pin that holds the blade in place. This had to be pushed out carefully. It will take some force as it is pressed in and will not fall out easily. Once this component is out, the old blade can be removed and the new one added in its place. Then the pin can be reinserted.

Click [HERE](#) for Tools for Chromatography Ordering Information

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)