



Remove a ferrule that is stuck in an HPLC system or column- How To

Sometimes a ferrule will become jammed inside a column or elsewhere in the HPLC system. When this happens, don't worry! The ferrule can be removed using our specially designed ferrule removal tool.

The tool works due to the direction of the threads, which go the opposite way from that of a standard drill bit. This allows the ferrule to be extracted out from the fitting.

To use, push against the ferrule with the tool while turning in a counter-clockwise direction until it is removed. If pushing is not enough, you may need to give it a light tap with a hammer to get it to bore into the ferrule.

Click [HERE](#) for Ferrule Puller Tool ordering information and pictures.

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com