



Are nylon membrane and syringe filters compatible with chloroform - FAQ

All nylon filters are not recommended for use with halogenated solvents such as chloroform or with strong acids (1N and stronger).

For these liquids, you should consider using PTFE filters.

Click [HERE](#) for ordering information and pictures of AQ brand of syringe filters.

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com