



Do the different tests in the HSQ Kit for HPLC use the same method conditions - FAQ

Yes, this is why it can be run overnight from an autosampler.

Simply set your conditions, program the system for performance qualification (PQ) validation and let it run. Use the following conditions:

Wavelength:	273nm
Injection Volume:	10µl (except for volume linearity)
Flow Rate:	1 mL/minute
Temperature:	Ambient

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com