



## How can I use the HSQ Kit to perform qualification for my specific analytes - FAQ

**The MicroSolv™ Qualification Kits are designed to qualify instruments and not analytes.**

What you may want be thinking of is system suitability which is different from performance qualification (PQ) and our kits are not for that use.

Once you qualify the instrument, it doesn't matter what types of analytes you are working with since the HSQ Kit™ qualifies the instrument performance itself. For example, once the instrument wavelength accuracy is qualified, you can have full confidence in this parameter regardless of what the UV spectra of your particular analytes look like. The full qualification of the various instrument parameters offered by this kit should exceed any regulatory agency's requirements for HPLC qualification.

Performance Qualification is a way to assure instrument accuracy and System Suitability assures precision of your method.

Click [HERE](#) for HSQ & PQ Kits ordering information and pictures.

Printed from the Chrom Resource Center  
Copyright 2025, All Rights Apply  
**MicroSolv Technology Corporation**  
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900  
Fax: (910) 769-9435  
Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)  
Website: [www.mtc-usa.com](http://www.mtc-usa.com)